

Floating Point Operation Error

Q: How do I solve the error “Floating Point By zero”?


A: Normally remove old document (un-install UM6-CAM software), and install again can work fine.

If still not working after re-install, it is possible because of the old document didn't remove completely. Please make sure NO **.ini** or **.txt** data left in UM6-CAM file.

Please follow below steps to remove software or file:

1. Open **UMX-CAM Viewer**. Go to **Tool > Open Data File > Enter” JPG File”**


(Data saving path ...\\MicroLinks Technology\\UM6-CAM\\...)

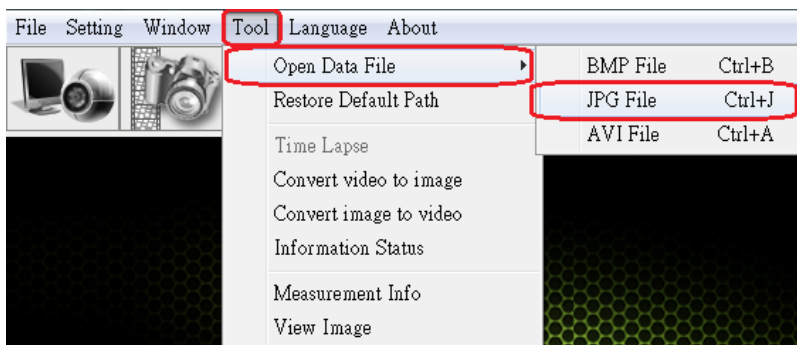

(Do not close the window)

2. Un-install UM6-CAM Software
3. Open the JPG folder and enter DATA folder. (Path: C:\\Users\\Administrator\\Microscope\\CAM6\\DATA)

Manually delete any **.ini** or **.txt** data in the folder. (ScaleData_Temp & UsersSysInfo)

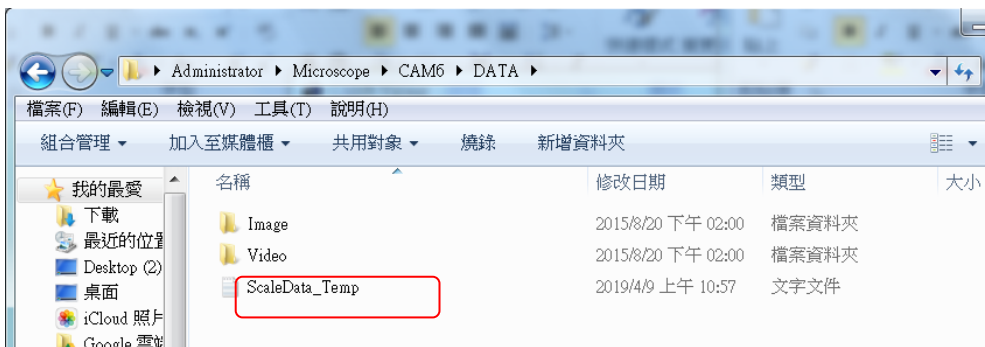

4. Install Software again.