



“UM Viewer” Software for IMB-06


- Setup the IMB-06 with ViTiny scope, and connect to computer.

Step 1. IMB-06 must initially boot up. Connect USB3.0 cable from PC/Laptop to  PC

Step 2. Connect mini USB cable from Microscope to  Device


Step 3. Connect mini HDMI cable from Microscope to  Device

Step 4. Connect HDMI cable from Monitor to  Monitor

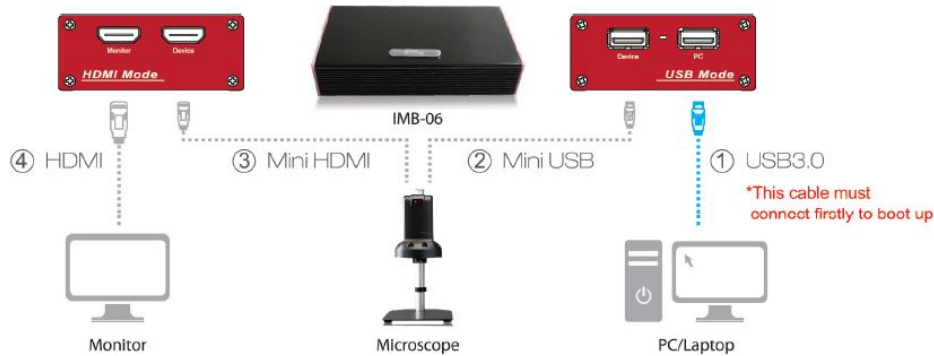


- Download “UM Viewer” software

1. Download link : http://www.vitiny.com/vitiny_en/ap_program.html
2. Open the UM Viewer and click “Connect” to start using software.

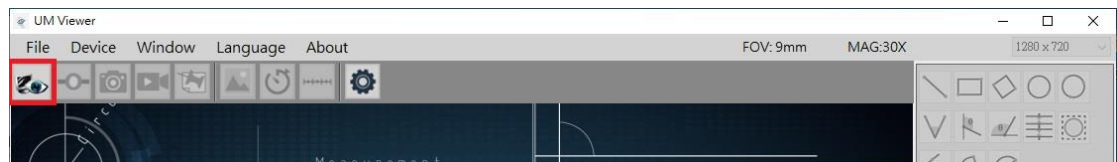


There are 4 advantage functions as below:


- When you press  on UM Viewer software. You can directly save photo in the PC and execute measurement etc. The remote controller become disable.




Control below functions like remote controller.



- **Auto calibration** : Press , please download video at <https://drive.google.com/file/d/1TByQy9Q3bcmt-asD-AUe11PcmzIVYJQI/view?usp=sharing>

- **Magnifier** :  measure helper.

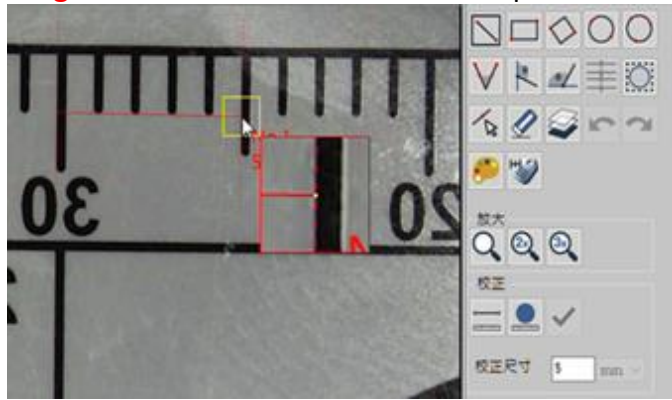

- **Edge detection** : 

