

MICROLINKS TECHNOLOGY CO.,LTD

User's Manual

Installing MICROSCOPE Device Operates Manual

2012/03/13 Update Ver.2.7A

Contents

1. Please insert CD in CD-Rom.....	3
2. Install/Uninstall MICROSCOPE application program	5
2.1 Install application program	5
2.1.1 Start setup program	5
2.1.2 Welcome to the MICROSCOPE Setup Wizard	5
2.1.3 Choose installation route.....	6
2.1.4 Now installing program	6
2.1.5 Install Completed	7
2.1.6 After installation completed.....	7
2.2 Uninstall MICROSCOPE AP	8
2.2.1 Choose MICROSCOPE AP's Uninstall as Fig. 2-8.....	8
2.2.2 Confirm to Uninstall MICROSCOPE.....	8
2.2.3 Uninstalling MICROSCOPE AP	8
2.2.4 Finish Uninstall.....	9
3. Uninstall MICROSCOPE Driver	10
3.1 Install driver	10
3.1.1 Start setup driver	10
3.1.2 Welcome to use USB video device setup wizard.....	11
3.1.3 Now installing.....	11
3.1.4 Install Completed	12
3.1.5 After installation completed.....	12
3.2 Uninstall MICROSCOPE Driver	13
3.2.1 Prepare to Uninstall. See Fig. 3-9	13
3.2.2 Confirm to remove the MICROSCOPE Driver	13
3.2.3 In process of uninstall	14
3.2.4 Finish Uninstall MICROSCOPE driver.....	14

1. Please insert CD in CD-Rom Slot

The package provides setup CD. If the setup CD is missing or damaged, please download setup files from <http://www.vitiny.com>. Please see the following steps to complete the MICROSCOPE installation.

- **Step 1: Please insert CD into CD-Rom.**
- **Step 2: Install MICROSCOPE application program.**
- **Step 3: Install MICROSCOPE driver.**

To uninstall MICROSCOPE AP and driver in your PC, please follow below 2 steps to complete the Uninstall.

- **Step 1: Uninstall MICROSCOPE AP.**
- **Step 2: Uninstall MICROSCOPE USB digital video driver.**

Please insert the product CD to the CD-Rom. The computer will read the CD automatically and will pop up a window shown below:



Fig.1-1 Setup window

If the shown setup window does not pop up, please open the files on the CD from your desktop and choose Autorun.exe file(see fig 1-2), then the same window (Fig 1-1) will pop up.



Fig.1-2 Autorun.exe icon

2. Install/Uninstall MICROSCOPE application program

Please select the Fig 1-1 Install program and setup application program button to install MICROSCOPE program.

2.1 Install application program

2.1.1 Start setup program

Click setup application program, after it will pop up (please select a language) window on Fig.2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click “Cancel” then click “OK”.

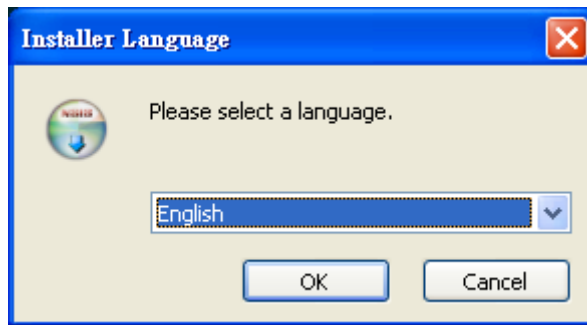


Fig.2-1 Setup language window

2.1.2 Welcome to the MICROSCOPE Setup Wizard

After choosing the language, go to welcome to the MICROSCOPE setup wizard window see Fig.2-2, then please click “Next” then continue. If want to cancel please click “Cancel” button. In the following figs, “name” represents MICROSCOPE’s application program name, “X.XXX.XXXx” represents the application program version.

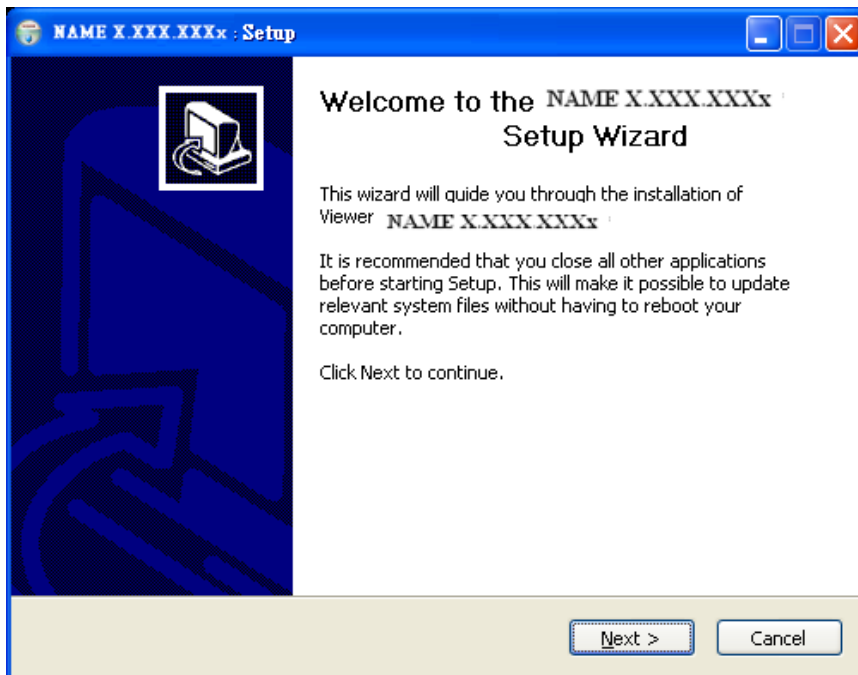


Fig. 2-2 Welcome to the MICROSCOPE setup wizard

2.1.3 Choose installation route

Next step please choose the installation route window see Fig.2-3. If want to change file route click “Browse” to change setup path, then please click “install” then start installing.

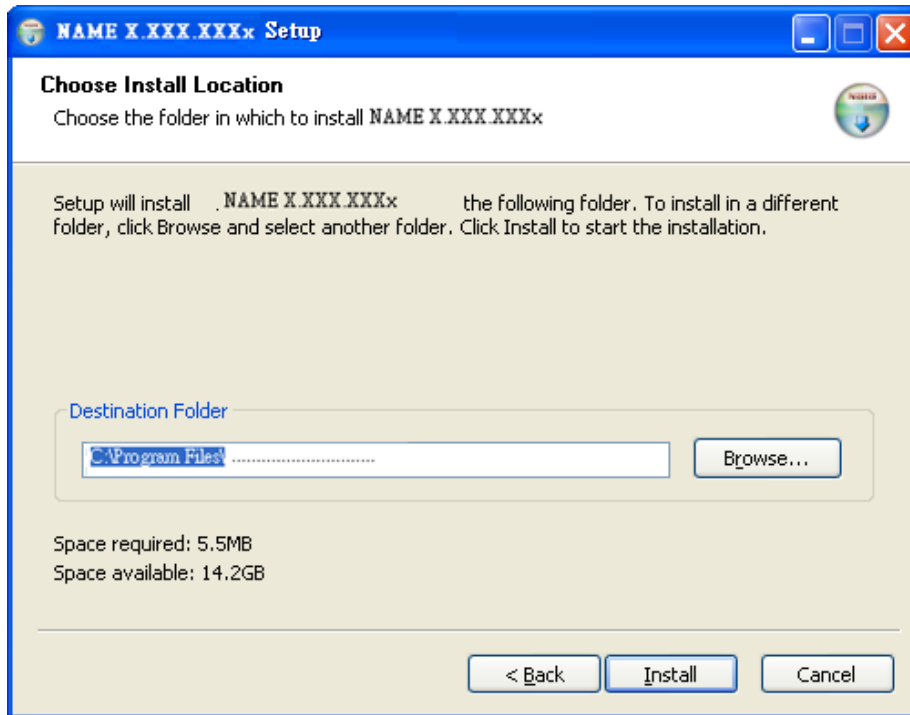


Fig. 2-3 Choose installation route

2.1.4 Now installing program

Continue install wizard will start installing program as directory Fig.2-4.

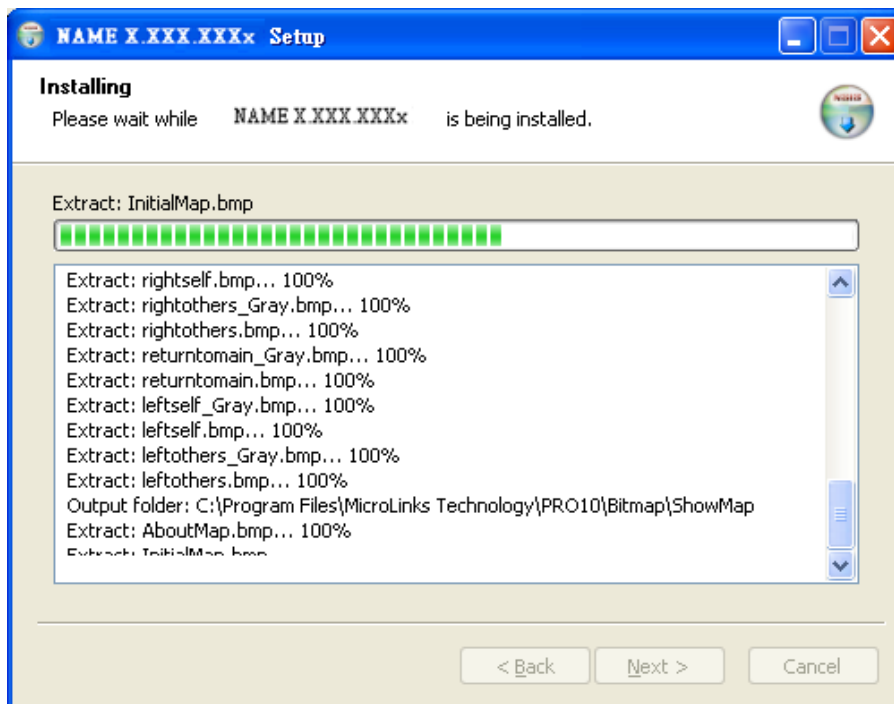


Fig.2-4 Now installing

2.1.5 Install Completed

When the MICROSCOPE program complete installation Fig.2-5. Please click “Finish”.

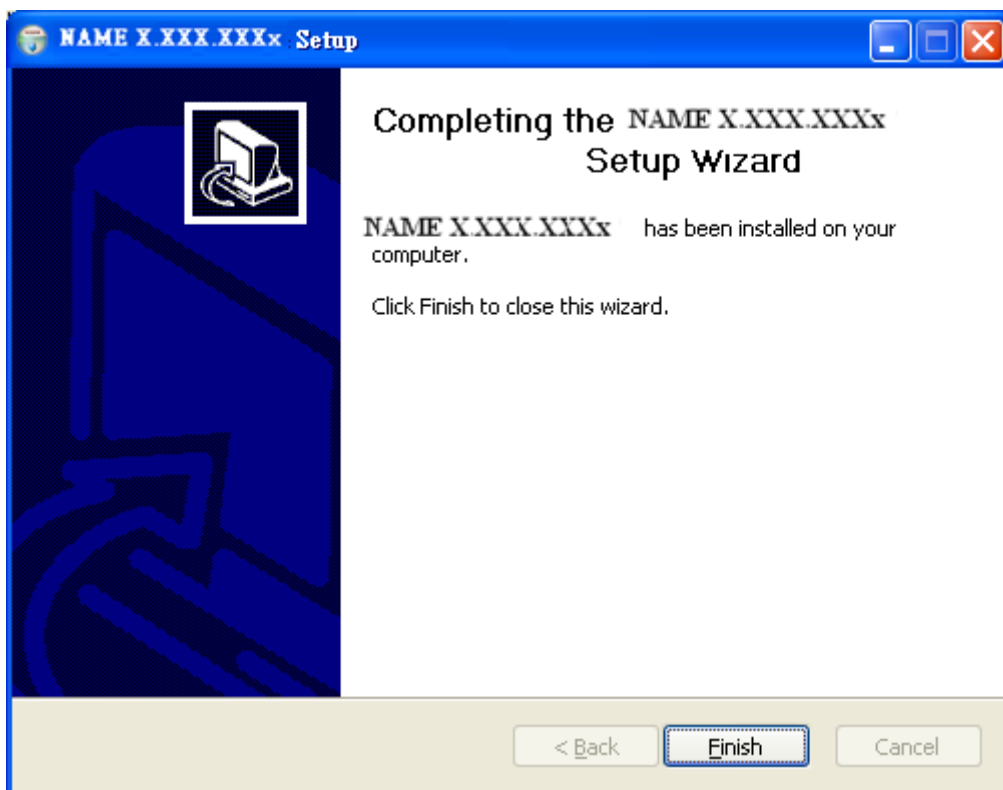


Fig.2-5 MICROSCOPE wizard install complete

2.1.6 After installation completed

MICROSCOPE application program shortcut will show on the Desktop as Fig.2-6.



Fig.2-6 MICROSCOPE shortcut

Or go “Start”-> All Programs-> MicroLinks Technology ->MICROSCOPE (see Fig 2-7).

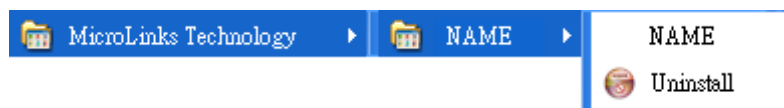


Fig.2-7 MICROSCOPE in All Programs

2.2 Uninstall MICROSCOPE AP

2.2.1 Choose MICROSCOPE AP's Uninstall as Fig. 2-8

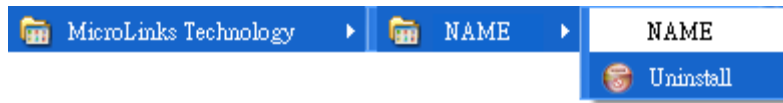


Fig.2-8 MICROSCOPE AP Uninstall

2.2.2 Confirm to Uninstall MICROSCOPE

Please make sure if you want to completely remove MICROSCOPE application program. See Fig.2-9

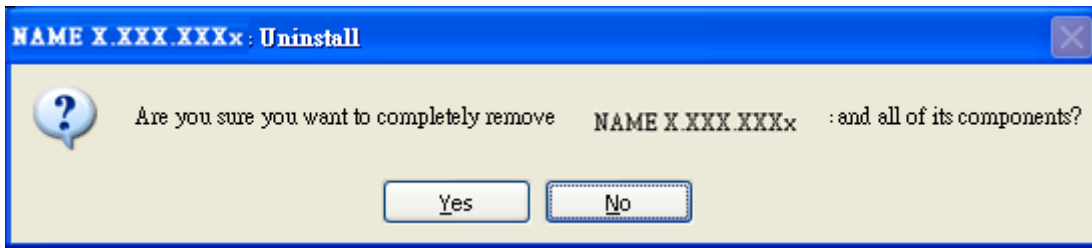


Fig.4-2

2.2.3 Uninstalling MICROSCOPE AP

In processing of uninstall the MICROSCOPE AP, see Fig.2-10

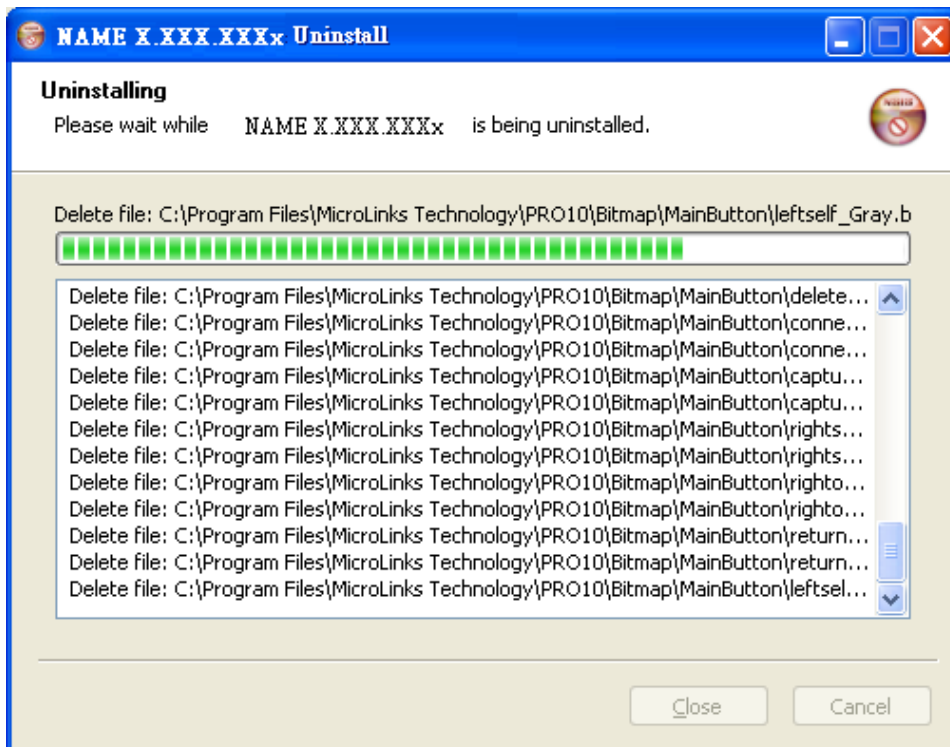


Fig.2-10 In process

2.2.4 Finish Uninstall

MICROSCOPE AP was successfully removed from your PC, please click OK to finish.



Fig. 2-11 Uninstall Completed

3. Uninstall MICROSCOPE Driver

The driver is optional one, if you install it, you can have more additional functions added to the MICROSCOPE AP.

Attention: Before install diver, please connect MICROSCOPE to the computer. If not, you can't install the program.

3.1 Install driver

3.1.1 Start setup driver

Please select Fig.1-1 to start Installing MICROSCOPE drive. Firstly please make sure to connect MICROSCOPE to a pc. Otherwise, it will pop up Fig.3-1. Note: Please use the USB cable from the gift box, do not use extension USB cable while connect microscope with a PC.

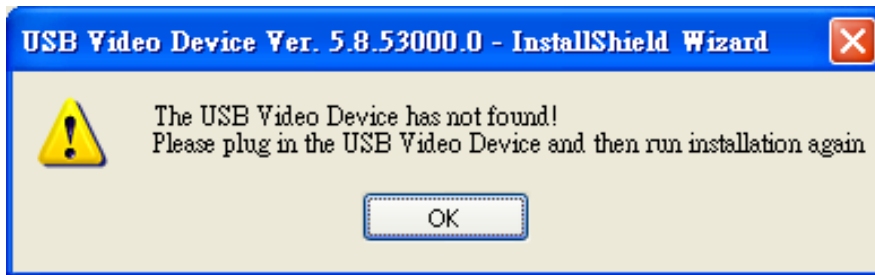


Fig.3-1 Not found MICROSCOPE step up

After click setup MICROSCOPE driver button, it will pop up Fig.3-2, now be ready to install the window. If cancel installation, please click "cancel" button.

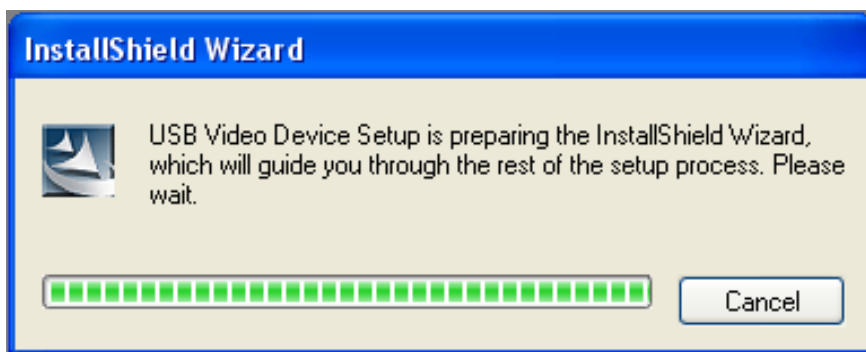


Fig.3-2 Getting ready to install

3.1.2 Welcome to use USB video device setup wizard

Ready to install, please follow up the steps.

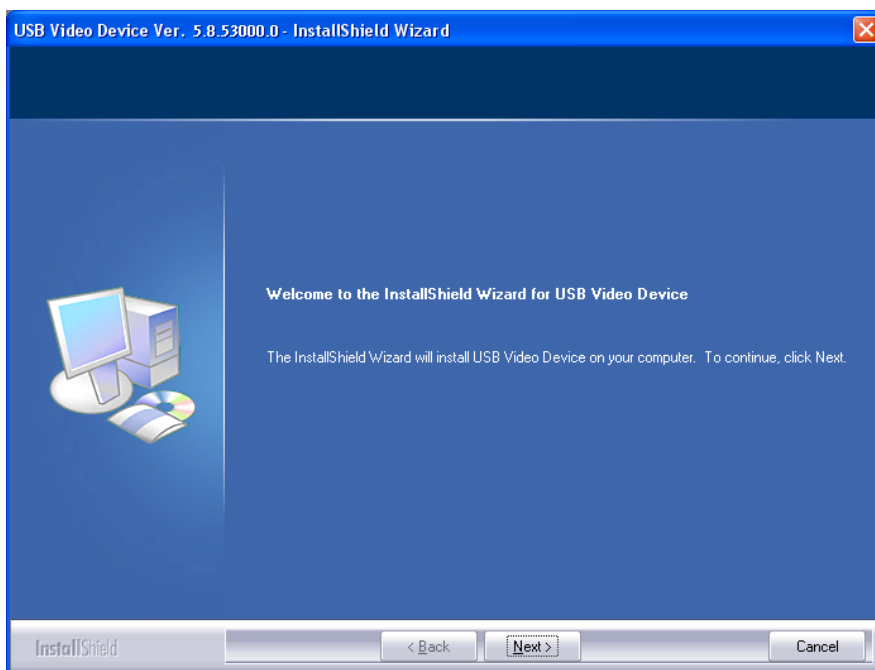


Fig.3-3 Welcome to use USB video device Install wizard.

3.1.3 Now installing

If you want to stop installation, please click "Cancel".

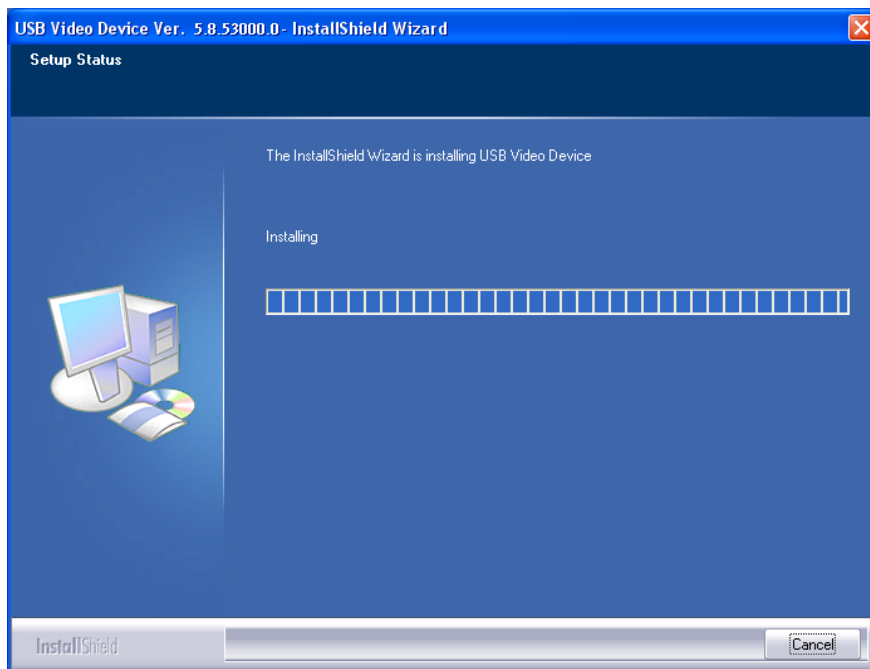


Fig.3-4 Now installing driver

3.1.4 Install Completed

Click “finish” to complete Driver Install program. See Fig 3-5.

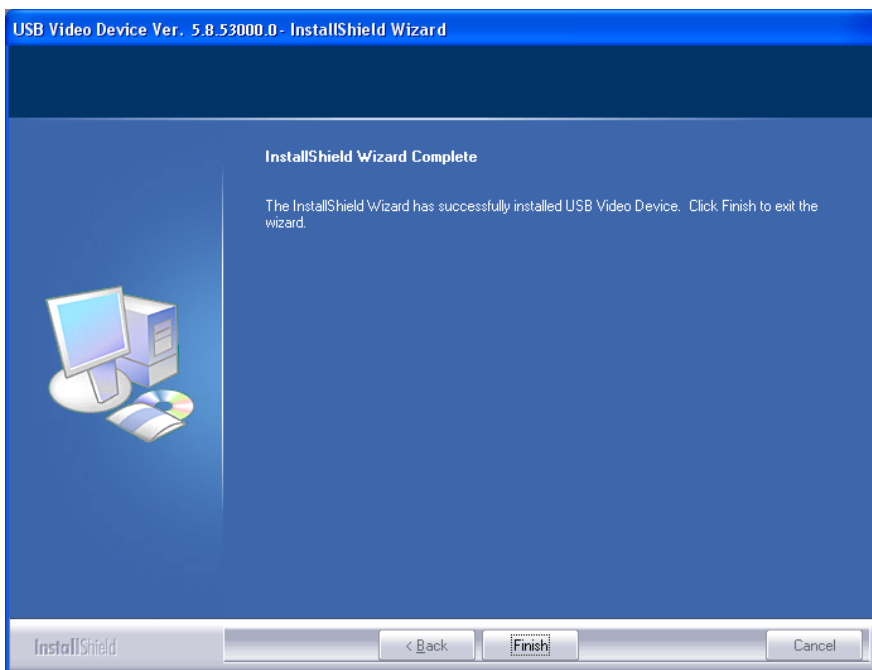


Fig.3-5 Driver installation completed

3.1.5 After installation completed

After installation is completed, the window on the right corner will show a video icon Fig.3-6. Start -> all programs will find Device Driver shown as Fig.3-7.



Fig.3-6 USB Video Device icon

MICROSCOPE can be executed from ICON AMCap as below Fig.3-6. However, AMCap is a simple tool for preview only and can't support image process as the MICROSCOPE AP.



Fig.3-7 USB Video Device is at all programs position.

3.2 Uninstall MICROSCOPE Driver

Choose MICROSCOPE driver icon to uninstall. See Fig.3-8



Fig. 3-8 Uninstall MICROSCOPE Driver

3.2.1 Prepare to Uninstall. See Fig. 3-9

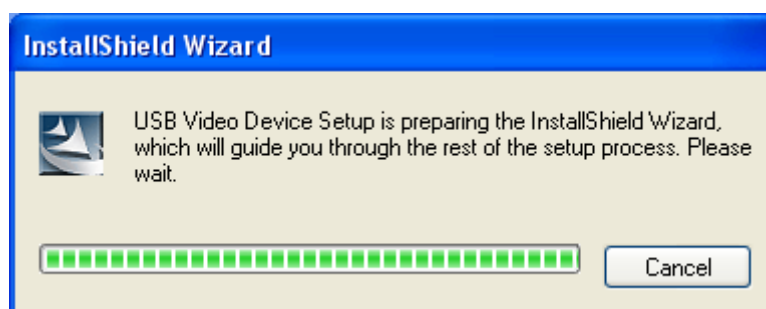


Fig. 3-9 Preparation

3.2.2 Confirm to remove the MICROSCOPE Driver

Please make sure if you want to completely remove the Driver. See Fig. 3-10

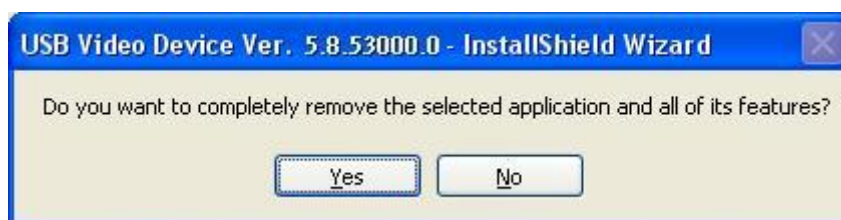


Fig. 3-10 Confirmation of uninstall

3.2.3 In process of uninstall

In processing of uninstall the Microscope Driver, see Fig.3-11

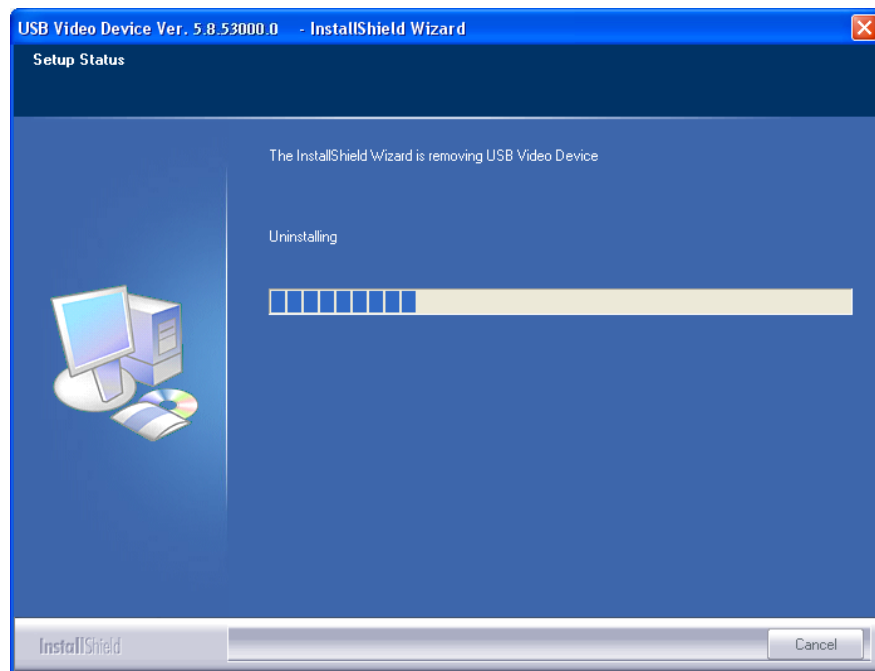


Fig.3-11 Uninstalling Driver

3.2.4 Finish Uninstall MICROSCOPE driver

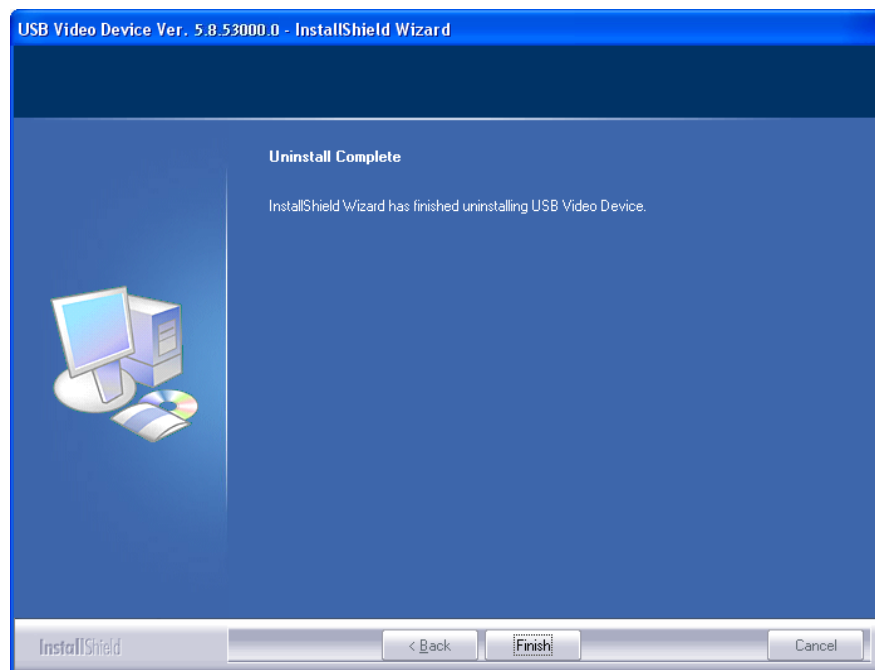


Fig.3-12 Finish uninstall