

MICROLINKS TECHNOLOGY CO.,LTD

# User's Manual

---

## Microscope Installation Manual

MLTC

2012/08/15 Update Ver.2.7A

## Content

<b>1. Please insert CD in CD-Rom.....</b>	<b>3</b>
<b>2. Install and uninstall Microscope Application Program.....</b>	<b>3</b>
2.1 Install application program .....	3
2.1.1 Start setup program .....	3
2.1.2 Welcome to the Microscope Setup Wizard .....	3
2.1.3 Choose installation route.....	4
2.1.4 Now installing program .....	5
2.1.5 Complete install .....	5
2.1.6 After installation completed .....	6
2.2. Install/ Uninstall Microscope AP .....	7
2.2.1 Choose MICROSCOPE AP's Uninstall as Fig. 2-8.....	7
2.2.2 Confirm to Uninstall MICROSCOPE.....	7
2.2.3 Uninstalling MICROSCOPE AP .....	8
2.2.4 Finish Uninstall.....	8

## 1. Please insert CD in CD-Rom

The package provides setup CD. If the setup CD is missing or damaged, please download setup files from <http://www.vitiny.com>. Please see the following steps to complete the Microscope installation.

- **Step 1: Please insert CD.**
- **Step 2: Install Microscope application program. Please refer to section 2.1.**

To uninstall Microscope AP in your PC, please follow below 2 steps to complete the Uninstall.

- **Step 3: Uninstall Microscope AP. Please refer to section 2.2.**

## 2. Install and uninstall Microscope Application Program

Please select the Fig.1-1 to start setup Microscope application program.

### 2.1 Install application program

#### 2.1.1 Start setup program

First, click “Install Microscope application program”, next, please select a language, see Fig. 2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click “Cancel” then click “OK”

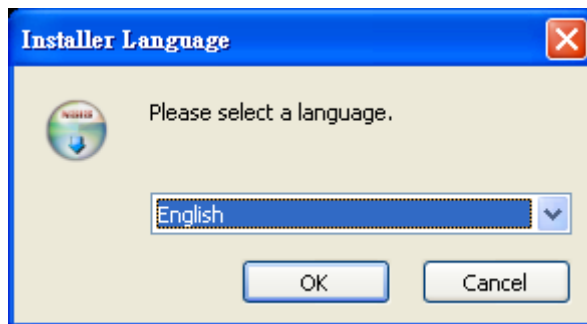


Fig 2-1seup language window

#### 2.1.2 Welcome to the Microscope Setup Wizard

After choosing the language, go to welcome to the MICROSCOPE setup wizard window see Fig.2-2, then please click “Next” then continue. If want to cancel please click “Cancel” button. In the folloing figs, “name” represents MICROSCOPE’s application program name, “X.XXX.XXXx” represents the application program version.

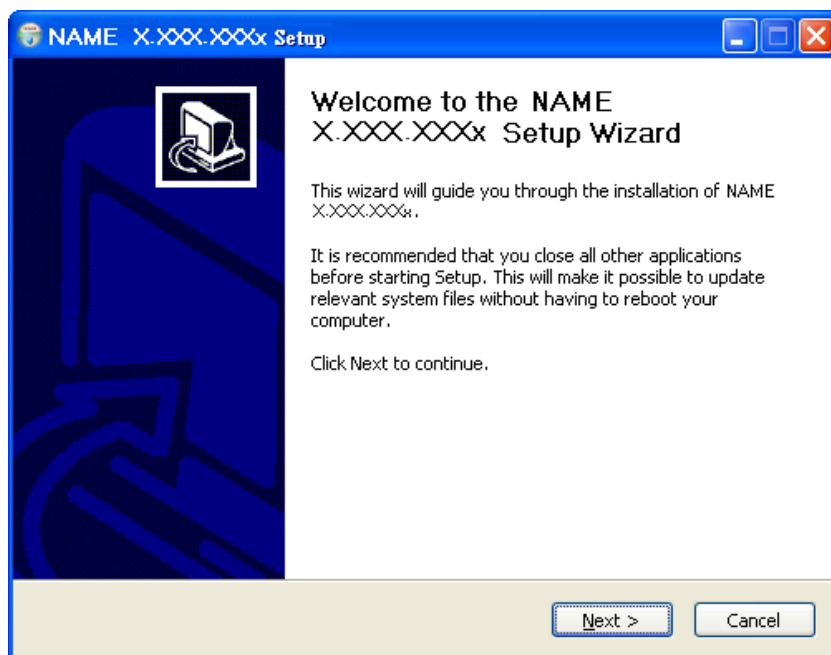


Fig 2-2 welcome to the Microscope setup wizard

### 2.1.3 Choose installation route

Next step please choose the installation route window see Fig.2-3. If want to change file route click “Browse” to change setup path, then please click “install” then start installing.

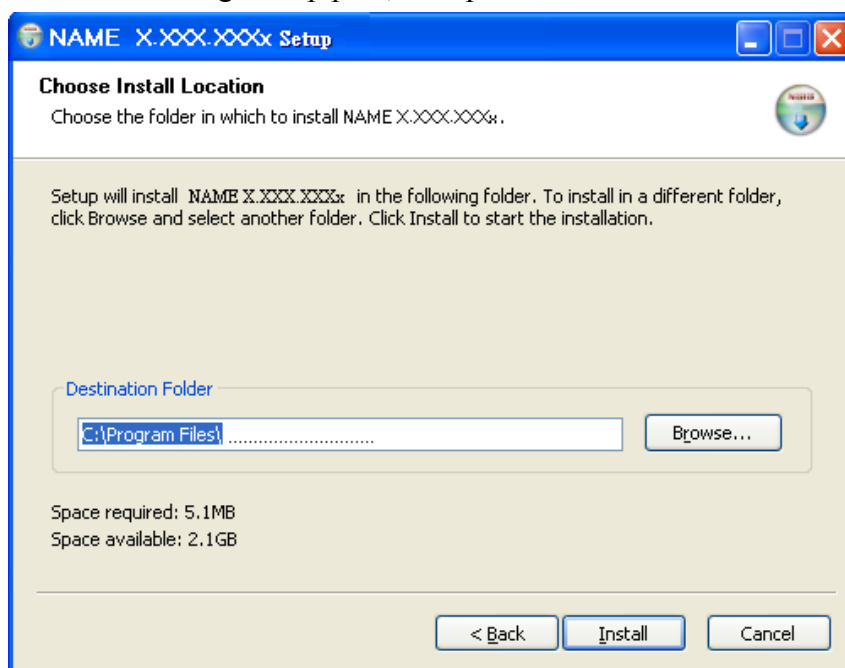


Fig 2-3 choose installation route

### 2.1.4 Now installing program

Continue install wizard will start installing program ad directory Fig 2-4.

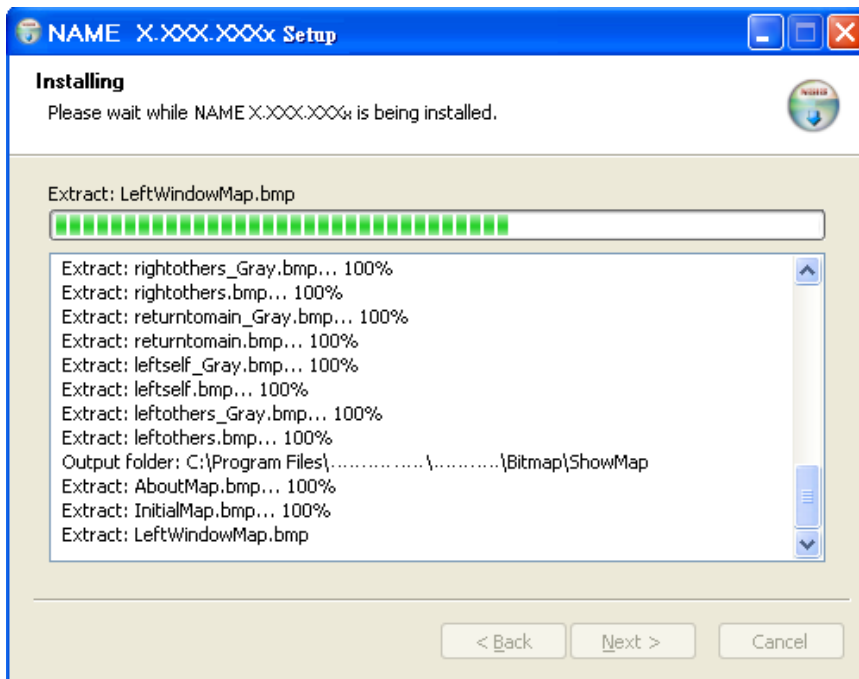


Fig 2-4 installing

### 2.1.5 Complete install

When the Microscope program complete installation (Fig 2-5), please click “Finish”.

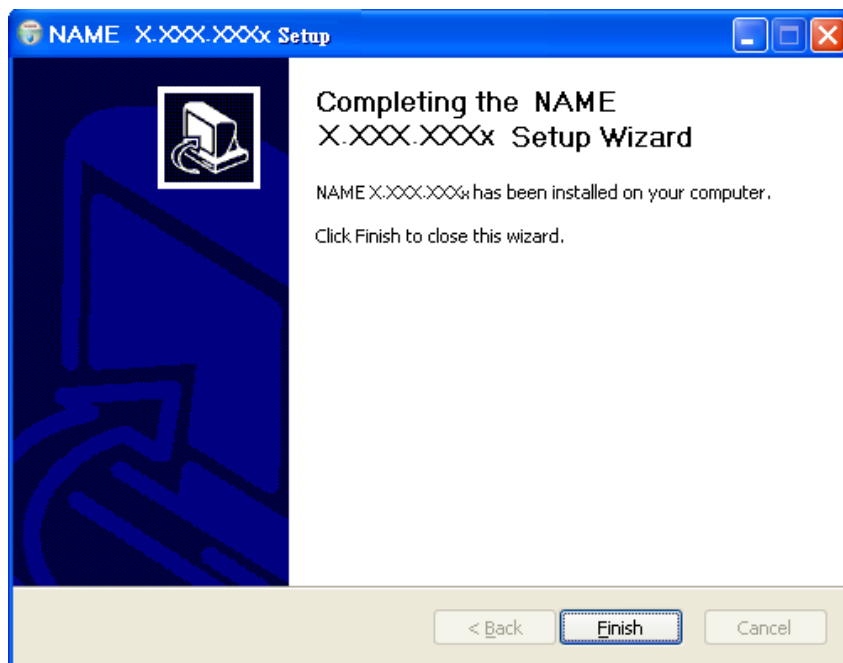


Fig 2-5 Microscope wizard install completed

### 2.1.6 After installation completed

MICROSCOPE application program shortcut will show on the Desktop as Fig.2-6.



Fig 2-6 Microscope shortcut icon

Microscope shortcut (Fig 2-6) will show on the Desktop. Or go “Start”-> All Programs-> MicroLinks Technology -> Microscope (see Fig 2-7).

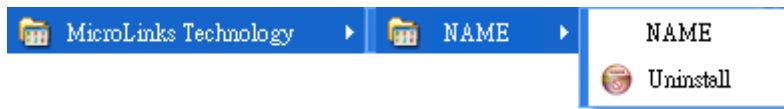


Fig 2-7 Microscope in All Programs

## 2.2. Install/ Uninstall Microscope AP

### 2.2.1 Choose MICROSCOPE AP's Uninstall as Fig. 2-8

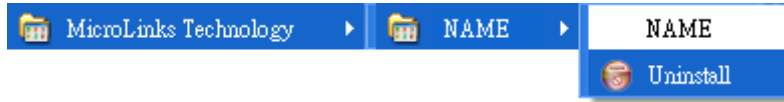


Fig.2-8 MICROSCOPE AP Uninstall

### 2.2.2 Confirm to Uninstall MICROSCOPE

Please make sure if you want to completely remove MICROSCOPE application program.  
See Fig.2-9

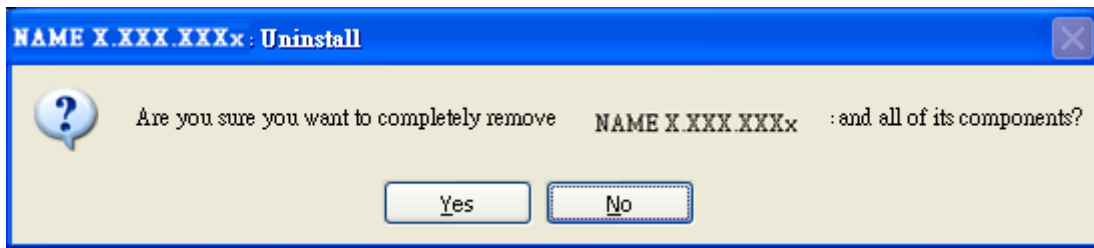


Fig.2-9

### 2.2.3 Uninstalling MICROSCOPE AP

In processing of uninstall the MICROSCOPE AP, see Fig.2-10

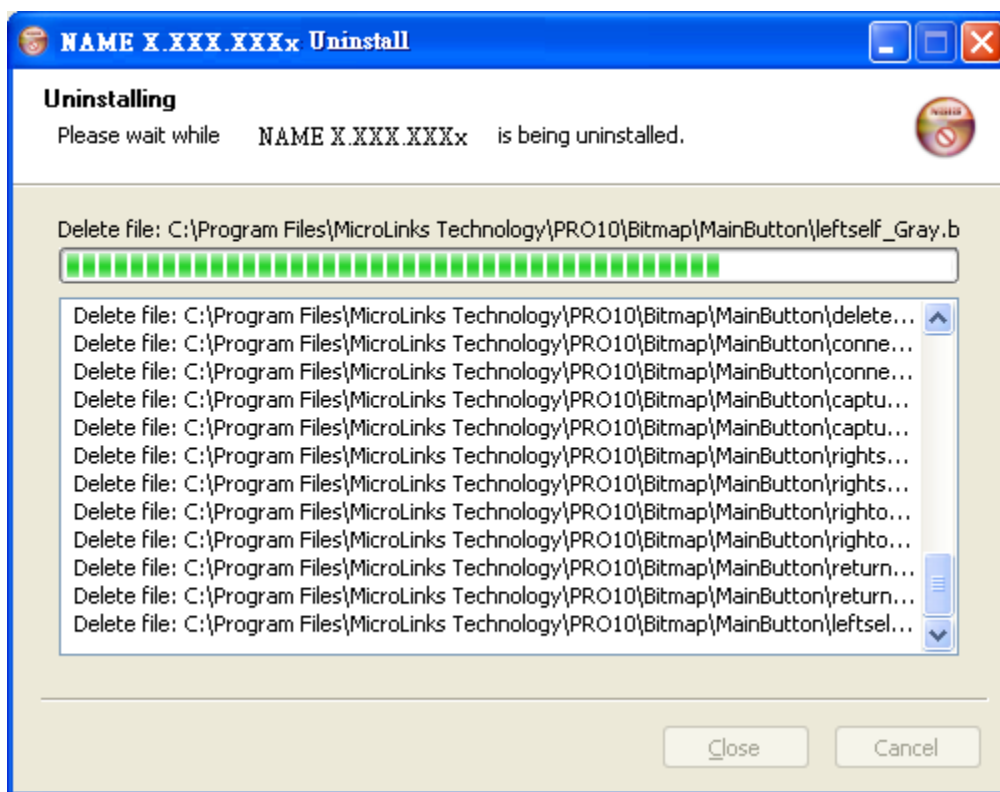


Fig.2-10 In process

### 2.2.4 Finish Uninstall

MICROSCOPE AP was successfully removed from your PC, please click OK to finish.

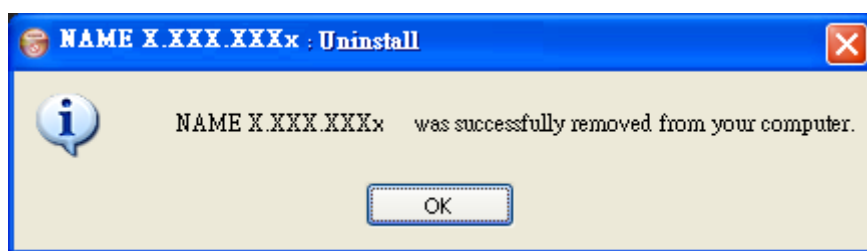


Fig. 2-11 Uninstall Completed