

MICROLINKS TECHNOLOGY CO.,LTD

# User's Manual

---

## Microscope Installation Manual

MLTC

2013/07/24 Update Ver.2.8A

## Content

1.	Introduction .....	3
2.	Microscope Application Program Installation and Uninstallation .....	3
2.1	Application Program Installation .....	3
2.1.1	Start the Setup Program .....	3
2.1.2	Welcome to the Microscope Setup Wizard .....	3
2.1.3	Choose Destination Folder .....	4
2.1.4	Installation Window .....	5
2.1.5	Complete Installation .....	5
2.1.6	Post Installation .....	6
2.2.	Application Program Uninstallation .....	7
2.2.1	Select the Uninstallation .....	7
2.2.2	Confirm to Uninstall Application Program .....	7
2.2.3	Uninstallation Window .....	8
2.2.4	Complete Uninstallation .....	8

# 1. Introduction

This guide outlines how to install and uninstall the microscope application program step-by-step. In the package, a setup CD should be provided. If the CD is missing or damaged, the setup files are available through download at <http://www.vitiny.com>.

## 2. Microscope Application Program Installation and Uninstallation

### 2.1 Application Program Installation

#### 2.1.1 Start the Setup Program

To begin, click “Install Microscope Application Program”, and then select a language from the drop-down menu as seen in Fig. 2-1. If you wish to cancel the setup please click “Cancel” and then click “OK”

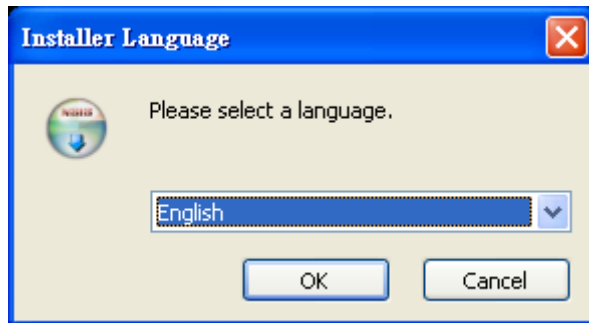


Fig 2-1 setup language window

#### 2.1.2 Welcome to the Microscope Setup Wizard

After selecting the language, the Microscope setup wizard will appear (see Fig.2-2). Click “Next” to continue or “Cancel” to exit. In the following figures, “name” represents the application program name and “X.XXX.XXXx” represents the application program version.

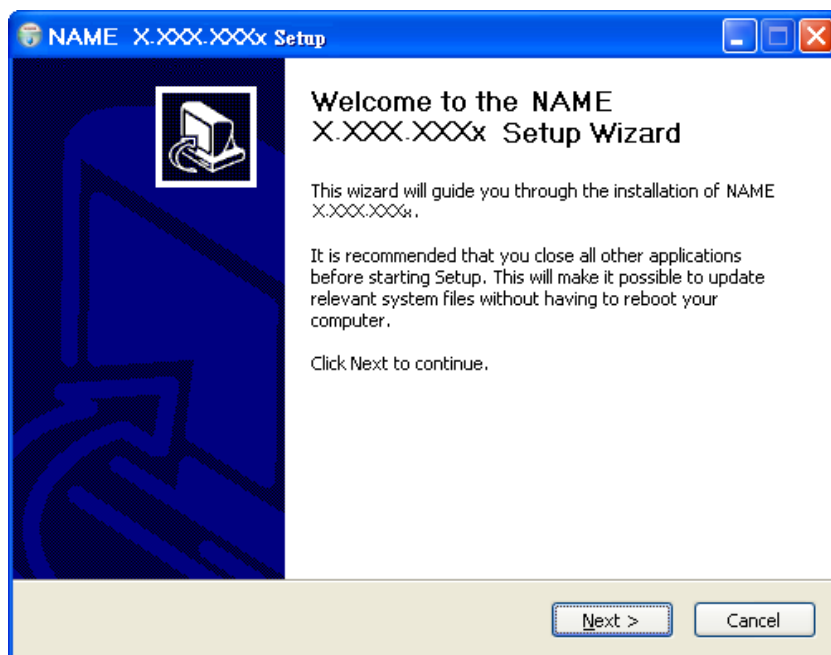


Fig 2-2 welcome to the microscope setup wizard

### 2.1.3 Choose Destination Folder

Choose the destination folder by clicking “Browse” and selecting a file destination (see Fig.2-3). Then click “Install” then start the installation.

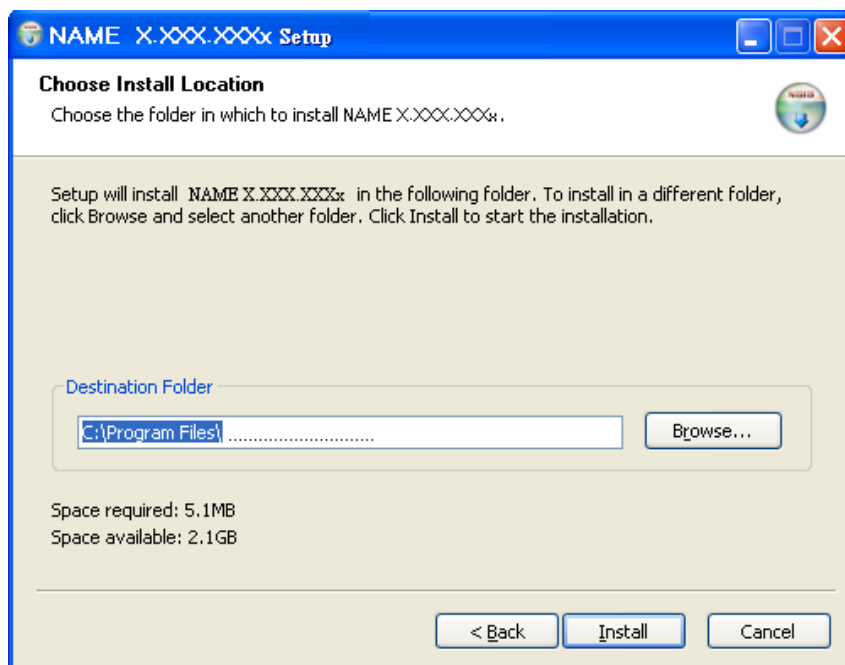


Fig 2-3 choose the destination folder

### 2.1.4 Installation Window

When the installation wizard (see Fig 2-4) is finished installing, click “Next”.

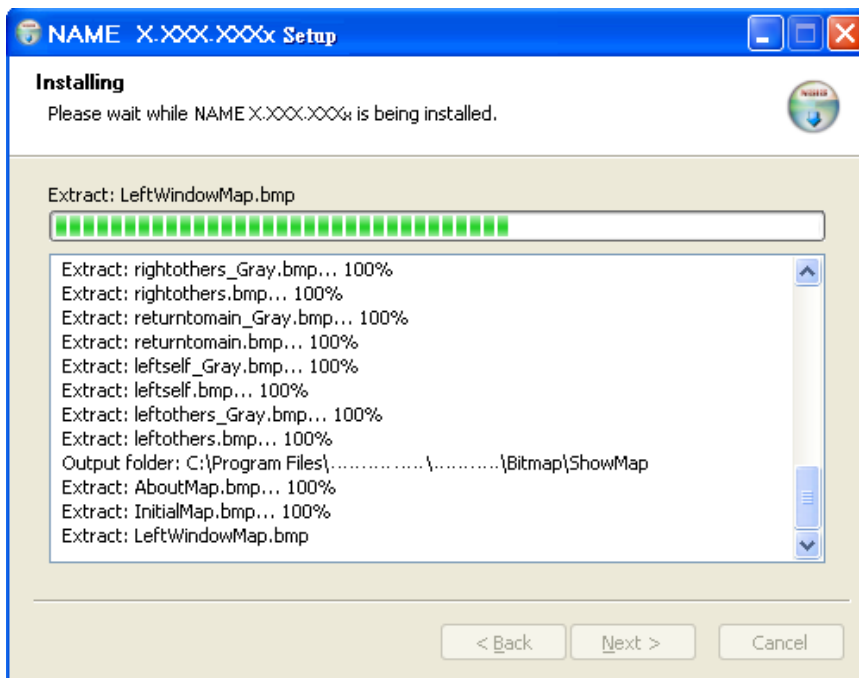


Fig 2-4 installing

### 2.1.5 Complete Installation

When the installation wizard has completed installation (Fig 2-5), click “Finish”.

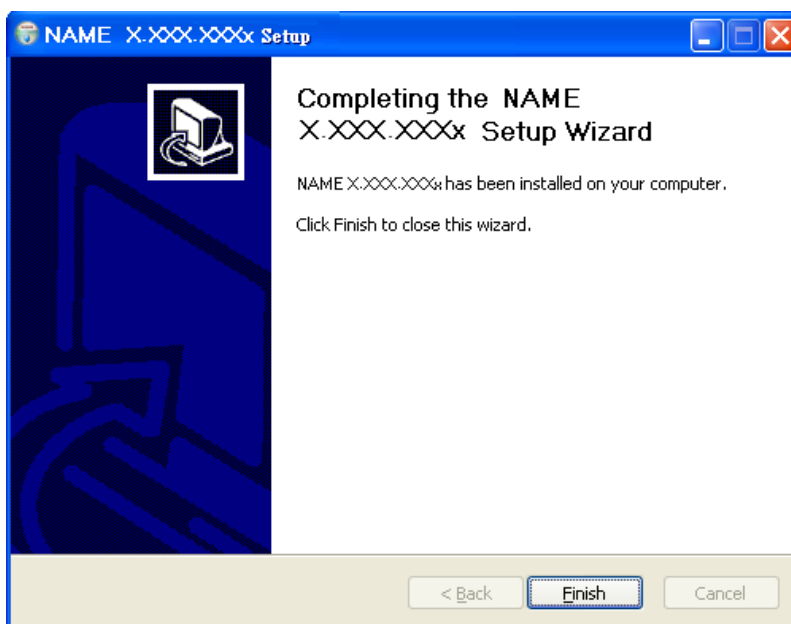


Fig 2-5 installation completed

### 2.1.6 Post Installation

The microscope application program is available through the desktop shortcut (Fig.2-6) or through “Start”-> All Programs-> MicroLinks Technology -> Microscope (Fig 2-7).



Fig 2-6 Microscope shortcut icon

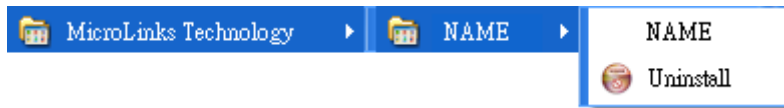


Fig 2-7 Microscope in All Programs

## 2.2. Application Program Uninstallation

### 2.2.1 Select the Uninstallation

Choose the microscope application “Uninstall” function as in Fig. 2-8

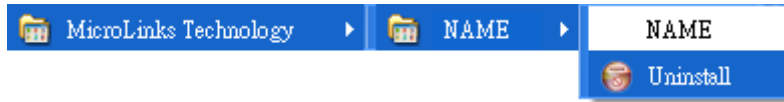


Fig.2-8 program uninstallation

### 2.2.2 Confirm to Uninstall Application Program

Select “Yes” if you want to completely remove the application program (Fig.2-9).

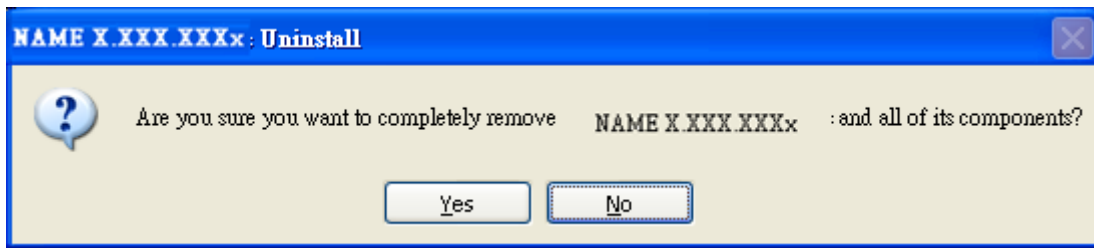


Fig.2-9 uninstallation confirmation

### 2.2.3 Uninstallation Window

When the installation wizard (see Fig 2-10) is finished installing, click “Close”.

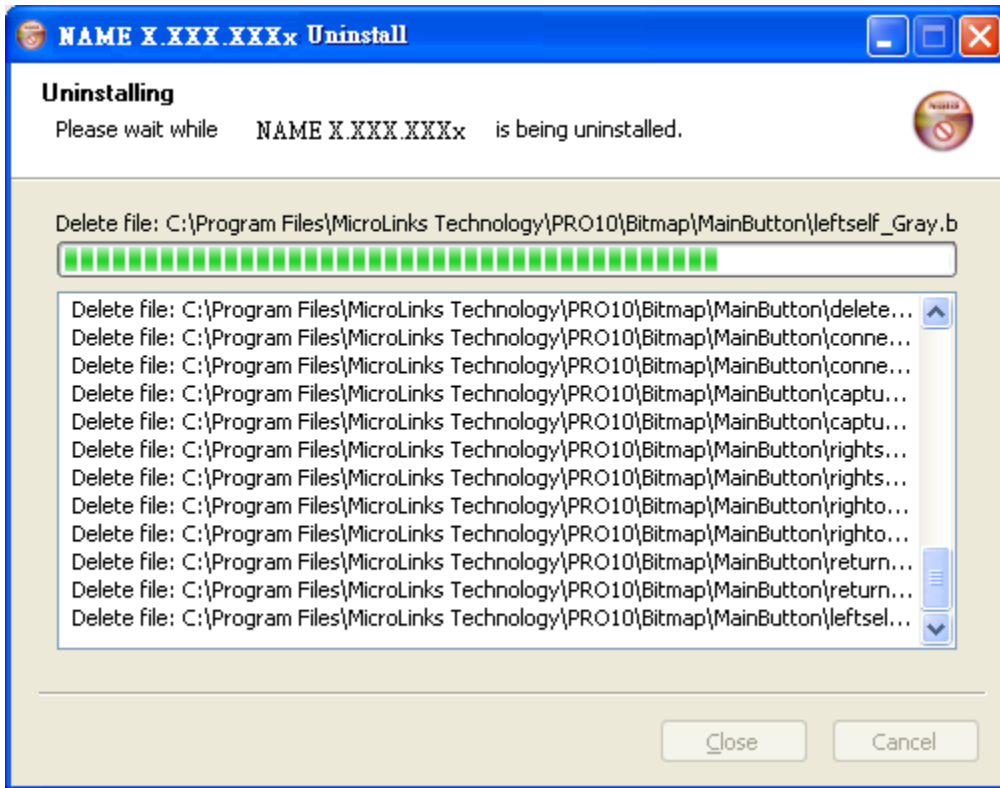


Fig.2-10 In process

### 2.2.4 Complete Uninstallation

Click “OK” to finish and exit.



Fig. 2-11 uninstallation completed